

Procedure for using UV Spectrophotometer (Lambda25)

1. Turn on the computer
2. Check the sample compartment to make sure that the beam paths are clear
3. Turn on the instrument. Do not open the compartment until the instrument is ready to use
4. Warm up for AT LEAST 10 minutes
5. Start UV Winlab (Click at the shortcut "Lambda25")
6. In the method window, click the tab "Wp"
7. Double click PKG825
8. Click "setup" to setup the instrument
9. Click "autozero"
10. Click "start"
11. Wait for the pop-up windows, Insert sample and click OK.
12. Repeat step 11 until all samples are measured.
13. Click "cancel"

Shutdown

1. Close down UV WinLab
2. Remove sample and reference
3. Clean the sample compartment and close it
4. Switch off the spectrophotometer
5. Turn off the computer